
**CKW BASE  
INSTRUCTION MANUAL**

**BASE CKW  
MANUAL DE INSTRUCCIONES**

**BASE CKW  
MANUEL D'INSTRUCTION**

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**BASE CKW  
MANUALE D'ISTRUZIONI**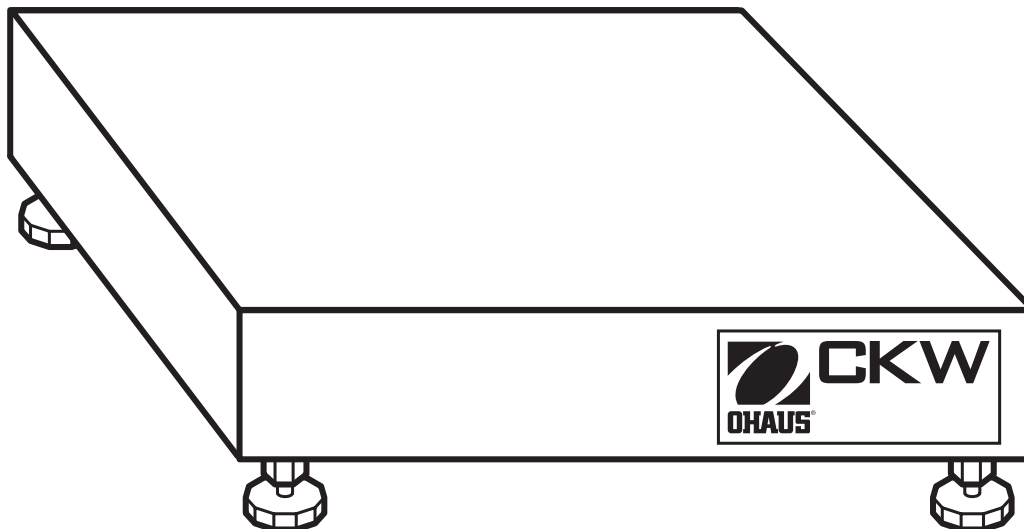

**Declaration of Conformity** We, Ohaus Corporation, declare under our sole responsibility, that the CE marked scale base models listed below are in conformity with the following directives and standards when used in combination with a CE marked indicator produced by Ohaus Corporation.

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Date: October 1, 2004

Ted Xia  
 President  
 Ohaus Corporation  
 Pine Brook, NJ USA

Urs Muller  
 General Manager  
 Ohaus Europe  
 Greifensee, Switzerland

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**ISO 9001 Registration**

In 1994, Ohaus Corporation, USA, was awarded a certificate of registration to ISO 9001 by Bureau Veritus Quality International (BVQI), confirming that the Ohaus quality management system is compliant with the ISO 9001 standard's requirements. On May 15, 2003, Ohaus Corporation, USA, was re-registered to the ISO 9001:2000 standard.

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# 1. INTRODUCTION

Thank you for purchasing a CKW Series product from Ohaus. Behind your purchase stands OHAUS, a leading manufacturer of precision Moisture Analyzers, Balances, Scales and Indicators. An Aftermarket Department with trained technicians is dedicated to provide you with the fastest service possible in the event your purchase requires servicing. OHAUS also has a Customer Service Department to answer any inquiries regarding applications and accessories.

This manual covers installation, operation and maintenance instructions for the Ohaus CKW Base. Please read this manual completely before installation and operation.

## 1.1. General Description

The Ohaus CKW Series is available as stand-alone scale bases or as complete weighing systems with the CKW-55 Indicator. The CKW Base is engineered for static checkweighing operations that also accommodates all types of weighing applications—from general purpose to washdown. With its easy-to-clean design to minimize food traps, food safety certification and rigid stainless steel construction, the CKW Base is ideally suited for years of use in harsh environments, and other wet-environment applications where sanitary requirements are high.

## 1.2 Features

- 4 Capacity Models: 3kg/6lb, 6kg/15lb, 15kg/30lb, 30kg/60lb
- 2 weighing pan sizes: 254 x 254mm (10" x 10") and 305 x 305mm (12" x 12")
- Resolutions: Maximum 1:6000 to 1:7500, Type-approved 1:3000
- Food Safety Certifications: NSF, USDA / AMS
- Supports HACCP compliance
- Stainless steel frame, pan and load cell (IP67 dust and water ingress protected)
- Adjustable non-slip rubber leveling feet with externally visible level indicator
- 6 rubber load pads and 6 down-stops for heavy-duty protection against shock and overload
- Manufactured under an ISO 9001 compliant quality system

## 1.3 Safety Precautions

For safe and dependable operation of this product, please comply with the following precautions:

- Operate the base only under ambient conditions specified in these instructions
- Ensure that the load cell cable does not pose an obstruction or tripping hazard
- Do not operate the base in hazardous environments or unstable locations
- Do not drop loads on the base
- Do not lift the base by the top plate; always lift from the bottom plate when moving the base
- Service should only be performed by authorized personnel

## 2. INSTALLATION

### 2.1 Unpacking

Unpack and inspect the product to make sure that all components have been included. The CKW Base package includes the following:

- CKW Base • Weighing Pan • Warranty Card • Instruction Manual

When purchased as a complete CKW scale, will also include:

- CKW-55 Indicator • Column Assembly

Check to ensure the product was not damaged during shipping. Immediately inform your OHAUS dealer in the unlikely event that you find damaged or missing parts.

Save the packaging material. Re-pack the product using original packaging materials to ensure safe storage or transport.

### 2.2 Assembly

#### 2.2.1 Weighing Pan

Place the weighing pan securely over the rubber load pads on the top plate of the base (first ensure that all load pads are seated fully in their respective mounting holes).

#### 2.2.2 Wiring Connections

Connect the load cell cable to an indicator using the wiring codes in Tables 2-1 and 2-2 below. Refer to the indicator manual for information about load cell connections, setup and calibration.

**Note:** When purchased as a CKW scale, the base is already pre-wired to the CKW-55 Indicator.

**TABLE 2-1. FOUR WIRE LOAD CELL CONNECTION  
(CKW3R, 6R and 15L)**

Function	Wire Color
+ Excitation	Green
- Excitation	Black
+ Signal	Red
- Signal	White

**TABLE 2-2. SIX WIRE LOAD CELL CONNECTION  
(CKW30L only)**


Function	Wire Color
+ Excitation	Green
- Excitation	Black
+ Signal	Red
- Signal	White
+ Sense	Blue
- Sense	Brown
Shield	Yellow

## 2.3 Selecting the Location

To ensure accuracy, proper performance and safety, locate and operate the base on a stable, level surface. Avoid locations with rapid temperature changes or excessive dust, air currents, vibrations, electromagnetic fields or heat.

Level the base by adjusting the four leveling feet until the bubble in the level indicator (located at the rear of the base) is centered. A wrench may be needed to loosen the locking nut above each leveling foot. When the base is level, retighten the locking nuts up against the base to lock each foot into place.

**Note:** Ensure that the base is level each time its location is changed.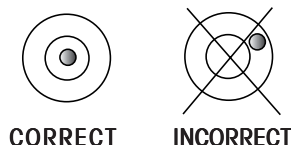

CORRECT

INCORRECT

## 3. MAINTENANCE

### 3.1 Cleaning

- The base components should be kept clean and free of excessive material build up.
- Avoid directly spraying the load cell with high-pressure water jets
- Water with a mild detergent may be used to clean the external and internal surfaces—do not use acids, alkalis, strong solvents or abrasive materials and chemicals
- The pan extends past the bottom plate to minimize entry of water or foreign material into the frame. As necessary, remove the pan to wipe or wash down the internal surfaces of the base.
- Removable components such as the weighing pan, leveling feet and rubber load pads (including column and CKW-55 Indicator when purchased as a scale system) may be removed and cleaned accordingly

### 3.2 Troubleshooting

Aside from installing components (Section 2.2) and leveling adjustments (Section 2.3), the CKW Base does not require any other adjustments as shipped from the factory.

Operational difficulties that may be encountered can often be traced to simple causes such as:

- Loose or incorrect wiring connections
- Obstructions to the weighing pan or base frame
- Unstable environments
- Incorrect calibration or setup of the indicator

If operational difficulties remain unresolved, contact your dealer or an authorized Ohaus Service Agent. Repairs should be made by authorized service personnel only.

For service assistance in the United States and Canada, you may call Ohaus Corporation toll free at (800) 526-0659 from 8 AM to 5 PM Eastern Time. An Ohaus Product Service Specialist will be available to help you.

## 4. TECHNICAL DATA

### 4.1 Drawings

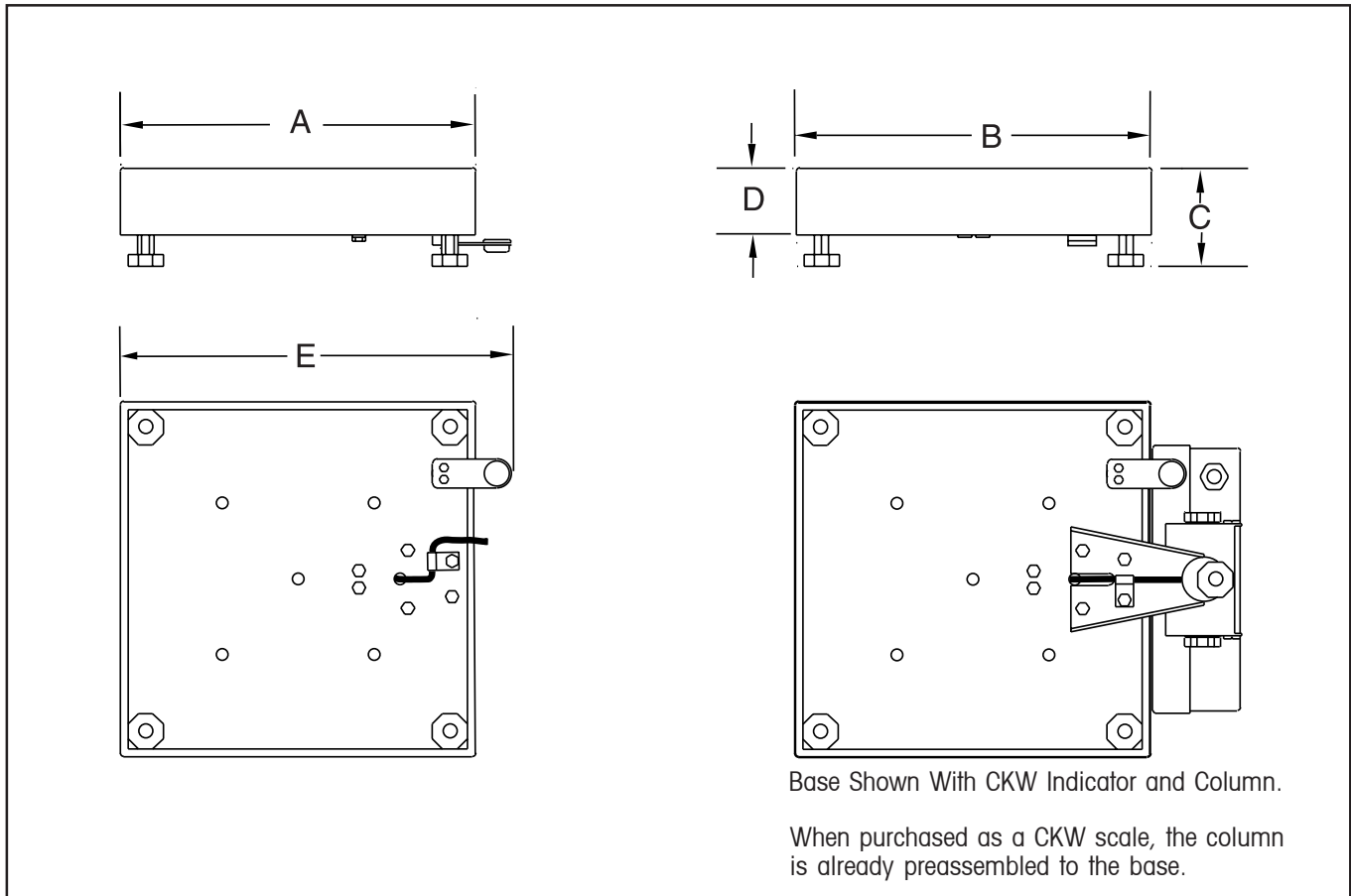
Figure 4-1. CKW Base Overall Dimensions.

### 4.2 Overall Dimensions

TABLE 4-1. DIMENSIONS.

Base	A	B	C	D	E
	Pan Depth	Pan Width	Height From Pan to Surface of Table*	Height of Pan	Depth Including Level Indicator
CKW3R, CKW6R	254 mm 10 in.	254 mm 10 in.	84 mm 3.3 in.	57.5 mm 2.26 in.	284.5 mm 11.20 in.
CKW15L	305 mm 12 in.	305 mm 12 in.	84 mm 3.3 in.	57.5 mm 2.26 in.	284.5 mm 11.20 in.
CKW30L	305 mm 12 in.	305 mm 12 in.	92 mm 3.6 in.	69.5 mm 2.74 in.	335.5 mm 13.21 in.

\* For leveling purposes, the feet may be extended up to an additional 11 mm (0.43 in.).

### 4.3 Technical Specifications

BASE	CKW3R	CKW6R	CKW15L	CKW30L
Capacity (max) x Readability (d)	3 kg x 0.0005 kg 6 lb x 0.001 lb	6 kg x 0.001 kg 15 lb x 0.002 lb	15 kg x 0.002 kg 30 lb x 0.005 lb	30 kg x 0.005 kg 60 lb x 0.01 lb
Maximum Resolution	1:6000	1:7500	1:7500	1:6000
Capacity (max) x Approved Readability (e)	3 kg x 0.001 kg 6 lb x 0.002 lb	6 kg x 0.002 kg 5 lb x 0.005 lb	15 kg x 0.005 kg 30 lb x 0.01 lb	30 kg x 0.01 kg 60 lb x 0.02 lb
Approved Resolution (n max)	1:3000	1:3000	1:3000	1:3000
Linearity	1d	1d	1d	1d
Repeatability (standard deviation)	1d	1d	1d	1d
Pan Dimensions (W x D x H)	254 x 254 x 57.5 mm 10" x 10" x 2.26 in.		305 x 305 x 57.5 mm 12" x 12" x 2.26 in.	305 x 305 x 69.5 mm 12" x 12" x 2.74 in.
Construction	304 Stainless Steel with Non-slip Rubber Leveling Feet			
Leveling Component	Externally Visible Level Indicator and Adjustable Leveling Feet			
Load Cell Capacity	6 kg	10 kg	30 kg	50 kg
Load Cell Cable	1m / 3ft. L x 4-wire	1m / 3ft. L x 4-wire	1m / 3ft. L x 4-wire	2m / 6ft. L x 6-wire
Load Cell Type	350 ohm, full bridge, Stainless steel			
Load Cell Excitation	5 to 15V DC/AC			
Load Cell Rated Output	2 mV/V			
Load Cell Protection	IP67			
Operating Temperature Range	-10°C to 40°C / 14°F to 104°F			
Relative Humidity	10% to 90%, non-condensing			
Height above sea level	Up to 4000 m			
Safe Overload Capacity	150% of rated capacity			
Product Weight	5.1 kg / 11.24 lb	6.2 kg / 13.67 lb	8.3 kg / 18.30 lb	9.3 kg / 20.50 lb
Shipping Weight	7 kg / 15.43 lb	8.1 kg / 17.86 lb	10.3 kg / 22.71 lb	11.3 kg / 24.91 lb
Food Safety Certifications	NSF, USDA / AMS			
Legal For Trade/Type Approval	<ul style="list-style-type: none"> <li>OIML: R60/2000-NL-02.14 / TC6046 r1 – 6, 10, 30 kg load cell</li> <li>R60/2000-NL1-03.24 / TC6120 r1 – 50 kg load cell</li> </ul>			

## 4.4 User-Replaceable Parts and Accessories

Part Number	Description
71170884	Pan, for CKW3R
71170885	Pan, for CKW6R
71170886	Pan, for CKW15L
71170887	Pan, for CKW30L
71170888	Rubber Load Pad, for Pan Support (Set of 6)
71170881	Leveling Foot (with locking nut)
71170126	Level Indicator Assembly (with mounting hardware)
71128010	Cable Clamp
71167972 <sup>1,3</sup>	Column Kit, for CKW-55, 229mm / 9" Fixed Height (All CKW Base models)
71154565 <sup>1,2,3</sup>	Column Kit, for CW-11, 350mm / 13.75" Fixed Height (All CKW Base models)
71154566 <sup>1,2,3</sup>	Column Kit, for CW-11, 680mm / 26.75" Fixed Height (CKW30L Base models only)
80251060 <sup>2,3</sup>	Column Adapter Kit, for connecting CW-11 column kits to CKW Bases

- Notes:**
- <sup>1</sup> All column kits listed above are of stainless steel construction. The CKW-55 Column Kit maintains the food use sanitary ratings.
  - <sup>2</sup> All CW-11 Column Kits require the Column Adapter Kit for proper connection to a CKW Base.
  - <sup>3</sup> If a column is later removed from the base, the original four bolts will need to be reinstalled at their original positions. This is required to maintain proper sanitary operation of the CKW base.

**LIMITED WARRANTY**

Ohaus products are warranted against defects in materials and workmanship from the date of delivery through the duration of the warranty period. During the warranty period Ohaus will repair, or, at its option, replace any component(s) that proves to be defective at no charge, provided that the product is returned, freight prepaid, to Ohaus.

This warranty does not apply if the product has been damaged by accident or misuse, exposed to radioactive or corrosive materials, has foreign material penetrating to the inside of the product, or as a result of service or modification by other than Ohaus. In lieu of a properly returned warranty registration card, the warranty period shall begin on the date of shipment to the authorized dealer. No other express or implied warranty is given by Ohaus Corporation. Ohaus Corporation shall not be liable for any consequential damages.

As warranty legislation differs from state to state and country to country, please contact Ohaus or your local Ohaus dealer for further details.

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